


The Confirmation Procedure of an upgrade period registered into HASP Key

1. Start HASP Info


You can confirm the upgrade period registered into the HASP key on the screen of Utility software called HASP Info. In order to invoke “HASP Info”, there are following two ways.

(a) Start from the Windows start

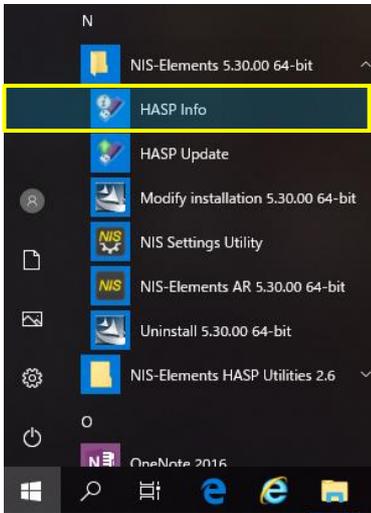

“HASP Info” of NIS-Elements is selected and started from the Windows start menu.

(b) Start from Help menu of NIS-Elements

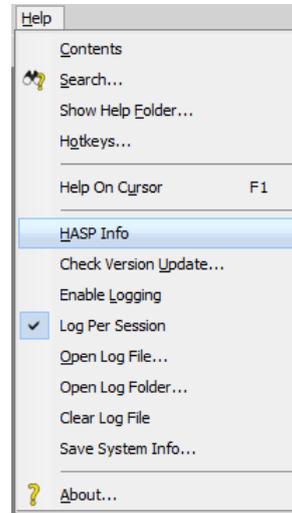

“HASP Info” is selected and started from Help menu of NIS-Elements after started NIS-Elements.

2. Confirmation of the upgrade period

The upgrade period displayed on a HASP Info screen is checked in the following procedures.

HASP SRM - Supported software and modules

HASP ID: 2710771B

Compatible with NIS-Elements 3.2 and higher

HASP 4 memory

		Version
Systems	NIS Elements AR	3.1
	NIS Elements BR	3.1
Modules	ND (6 dimensions)	

HASP SRM memory

		Version
Systems	NIS Elements AR	3.2
	NIS Elements BR	3.2
Modules	ND (6 dimensions)	

Activation Date: August 02, 2010 (1) Activation Date

Upgrades available until September 02, 2011 (2) Upgrade Period

The following two information is shown at the bottom of “HASP Info”.

- (1) Activation Date
- (2) Upgrade Period

If the release date of each version of NIS-Elements is before the date which is displayed in the upgrade period, you can upgrade to the version.

If the release date is after the date which is displayed in the upgrade period, to upgrade the version, you need to purchase the SUA option for a required period, and need to extend an upgrade period.*1

*1: If you have NIS-Elements D and the activation date was set before the delivery date, please contact your place of purchase for assistance. There is a possibility that your product was shipped after the free upgrade period had expired. By applying the SUA free of charge, we will ensure that the planned release date of Ver7.0 is within the SUA validity period. This will allow you to upgrade to Ver7.0 at no additional cost.

3. Confirmation for release date of NIS-Elements

You can check "release date" of each version of NIS-Elements at download for NIS-Elements.
In the example of the following figure, the date is displayed in red box is release date of NIS-Elements.

NIS-Elements AR

Imaging Software NIS-Elements AR Ver5.00.00 for 64bit edition

Login / New Registration
To download software and/or firmware, user registration is required.

> [User Registration](#)

Imaging Software **NIS-Elements AR Ver5.00.00 for 64bit edition**

Use Condition

1. Conditions of Use
 - This upgrade can be applied to the following version:
NIS-Elements AR Ver5.00.00 64bit.
 - Check "HASP Info" to confirm below:
"Upgrades available until May 24, 2017" or later.
 - [For details, please refer to "Confirmation Procedure of an upgrade period". \(PDF:96KB\)](#)
2. Installation

Release date:26/05/2017

Summary

The following figures are the summaries about the upgrade period which was registered into the HASP Key, Release date, and purchase of SUA option.

