

NIS-Elements L

License Manual

1st Release

Contents

1	Introduction	3
2	Preparing.....	3
3	The procedure	3
3.1	Online transfer (in the case that the both PCs are connected to the Internet)	3
3.2	Offline Transfer (in the case that the both PCs are not connected to the Internet)	5
4	[Important] Caution	8

1 Introduction

The license transfer is necessary to change the PC where NIS-Elements L is operated. This document describes the procedure.

2 Preparing

Please prepare the following items for license transfer.

- PC where NIS-Elements L activation has been done
- PC to which NIS-Elements L is to be used after this.
- Internet environment.


If the PCs are not connected to the Internet, please prepare another PC connected to the Internet.

3 The procedure


3.1 Online transfer (in the case that the both PCs are connected to the Internet)

- 1) Evacuating the license (Please operate this at the PC from which license is transferred)


Please start licmgr.exe, which exists the NIS-Elements L install folder (Default: C:\Program Files\NIS-Elements L). The following window is displayed.

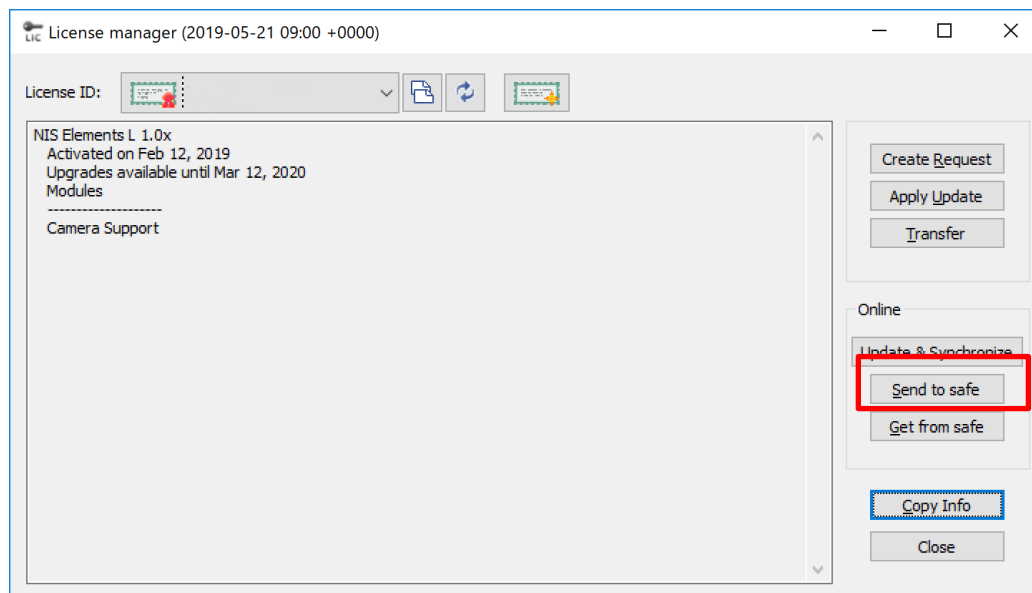

Please click "Save to safe" button. Then, input any name and password and click "Next" button. By this operation, the license is evacuated to the license server.

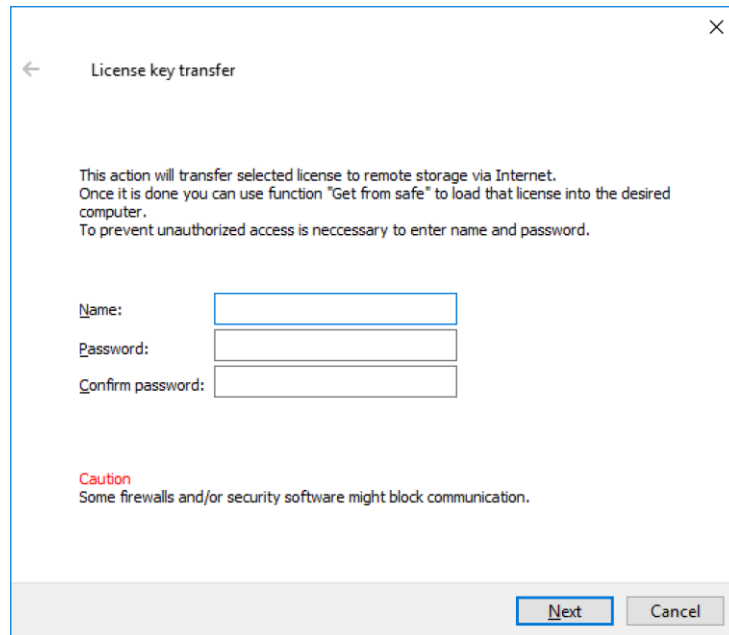


- 2) Loading the license (Please operate this at the PC to which the license is transferred)
Please start licmgr.exe, which exists the NIS-Elements L install folder (Default: C:\Program Files\NIS-Elements L).

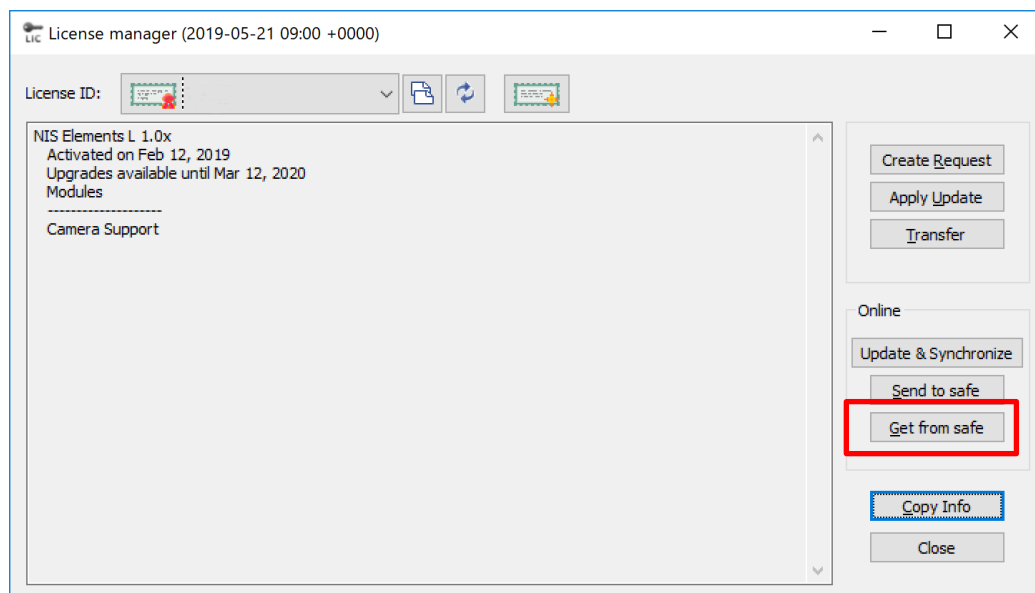

Please click "Get from safe" button. Then, input the name and password which are set in the previous step and click "Next".

After this, the license transfer is completed.

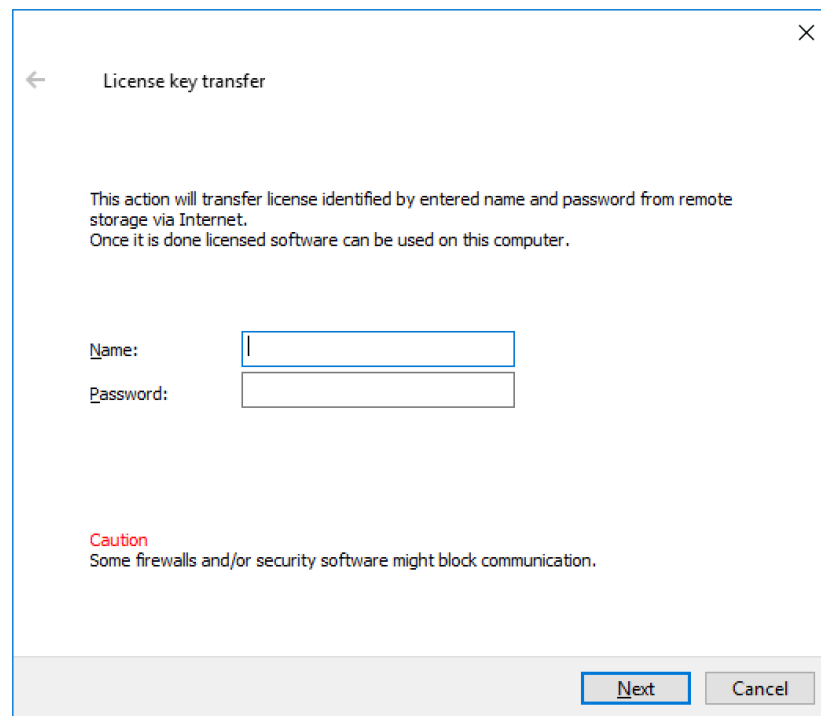

A screenshot of a Windows dialog box titled "License key transfer". The dialog has a standard Windows window border with a close button (X) in the top right corner. Inside the dialog, there is a back arrow button in the top left. The main text reads: "This action will transfer license identified by entered name and password from remote storage via Internet. Once it is done licensed software can be used on this computer." Below this text are two input fields: "Name:" and "Password:". At the bottom right, there are two buttons: "Next" and "Cancel". A "Caution" message in red text is located above the buttons, stating: "Some firewalls and/or security software might block communication."


3.2 Offline Transfer (in the case that the both PCs are not connected to the Internet)

In this way, another PC which is connected to the Internet is needed.

- 1) Creating the fingerprint of new PC (Please operate this at the PC to which the license is transferred)

Please start licmgr.exe, which exists the NIS-Elements L install folder (Default: C:\Program Files\NIS-Elements L). The following window is displayed.


Click "Create Request" button and save the c2l file. Then, move that c2l file to the PC from which the license is transferred.

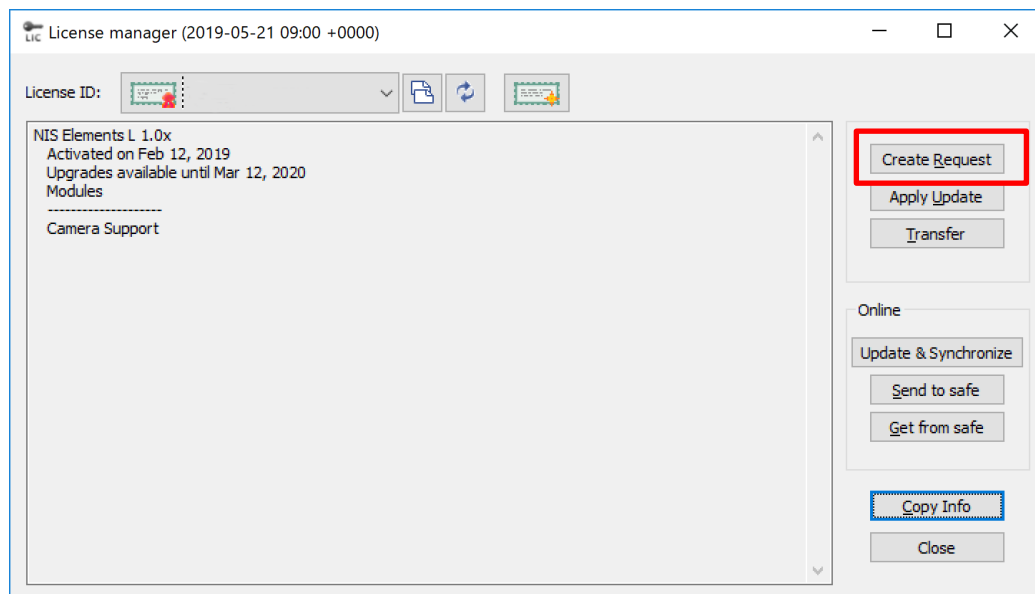

- 2) Creating the intermediate transfer file (Please operate this at the PC from which the license is transferred)

Please start licmgr.exe. Then, click "Transfer" button and have the c2l file created at the previous step loaded.

Then, the intermediate transfer file (h2l file) is created. Save it and move it to the PC which is connected to the Internet.

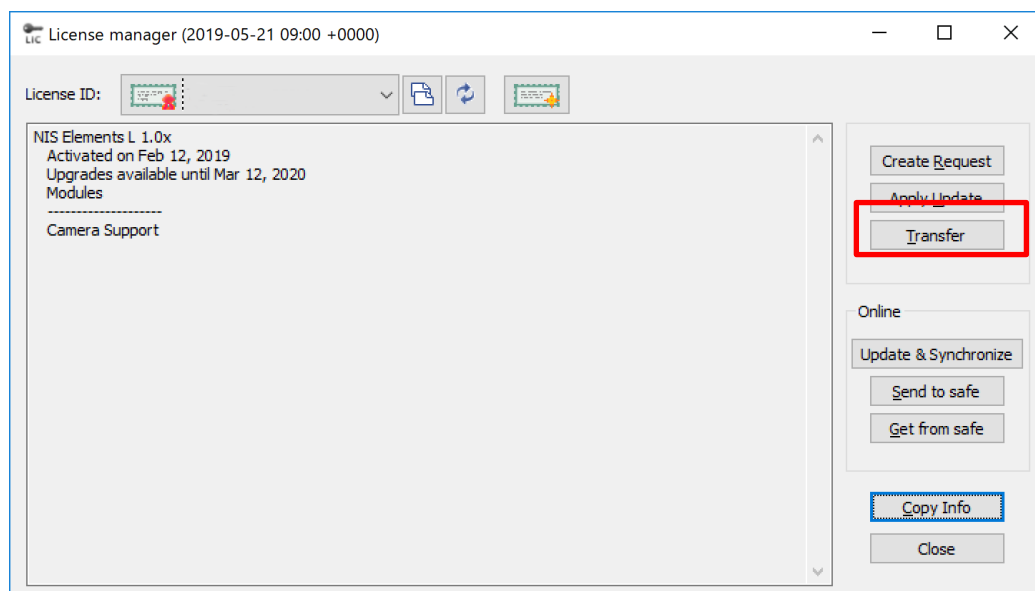


- 3) Creating the transfer file. (Please operate this at the PC which is connected to the Internet)

Please access to the NIS-Elements L Transfer site using a browser.

NIS-Elements L Transfer site :

<https://www.nis-activation.com/transfer/>

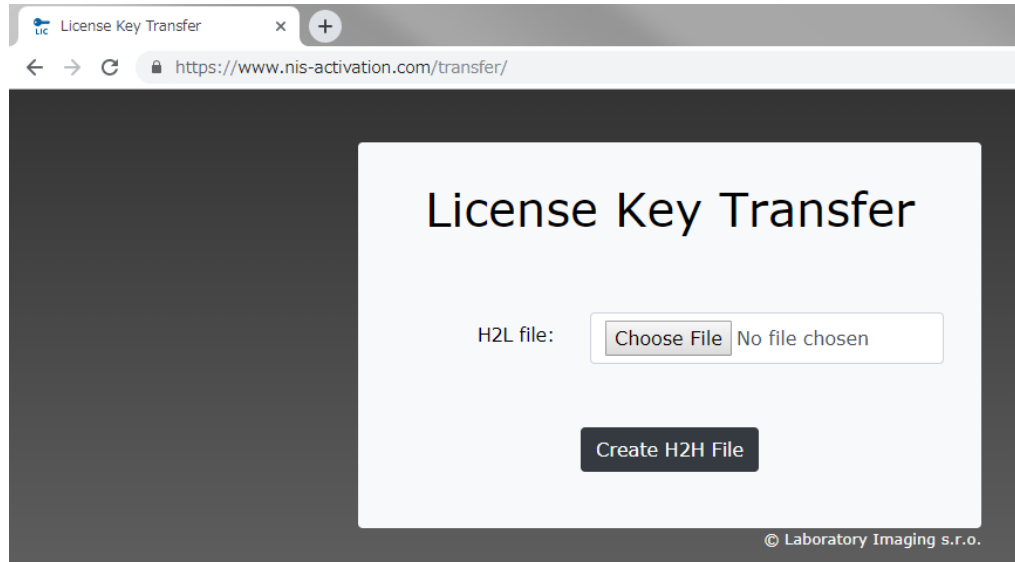


Please click the "Choose File" button and have the h2l file created at the previous step loaded.

Then, the transfer file (h2h file) is created. Please save it and move it to the PC to which the license is transferred.

- 4) Applying the transfer file (Please operate this at the PC to which the license is transferred)

Please start licmgr again, click "Apply Update" button and have the h2h file created at the previous step loaded.

Then, license transfer is completed.

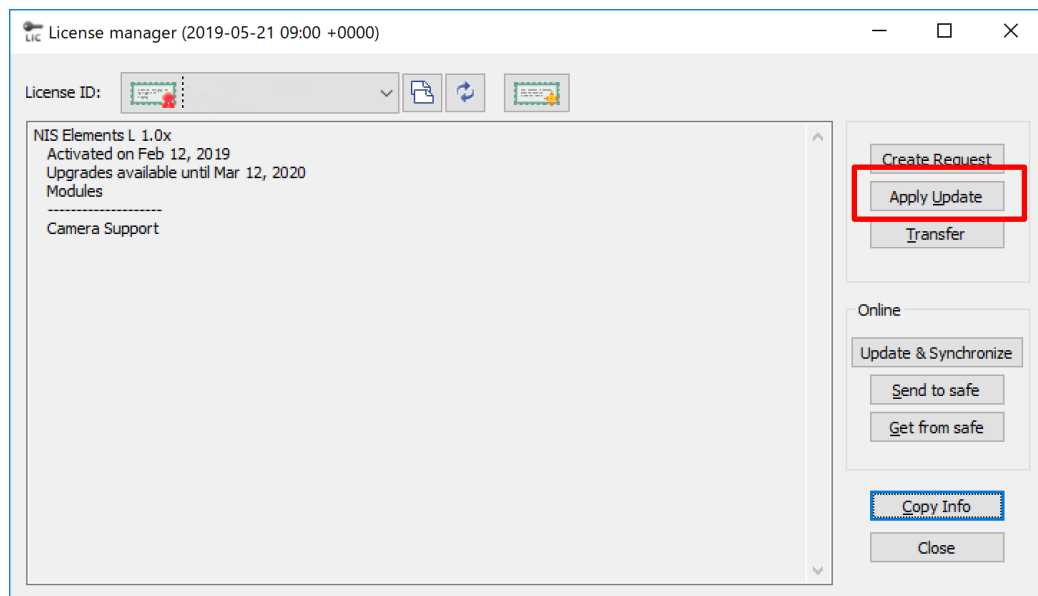

4 [Important] Caution

When you want to initialize the PC where license activation is done, please follow the below procedure.

1. Have the PC connected to the Internet.
2. Evacuate the license (see chapter 3.1 step1)
3. Initialize the PC
4. Have the PC connected to the Internet again
5. Load the license (see chapter 3.1 step2)

Without this procedure, the license is expired and NIS-Elements L cannot be used by that license after this.

End