

NIS-Elements L

Activation Manual

2nd Release

Contents

1	Introduction	3
2	Preparing.....	3
3	The procedure in the case that the PC where NIS-Elements L is installed is connected to the Internet.....	3
3.1	Online activation	3
3.2	In the case that online activation is failed	4
4	In the case that the PC where NIS-Elements L is installed is not connected to the Internet	6
4.1	Creating the request file (C2L)	7
4.2	Creating the activation file (V2L)	7
4.3	Activating the license.....	8
5	[Important] Caution	8

1 Introduction

The license activation is necessary to start NIS-Elements L. This document states the procedure of license activation.

2 Preparing

Please prepare the following items before activation.

- PC where NIS-Elements L has been installed
- License card
- Internet environment


If the PC where NIS-Elements L is installed cannot be connected to the internet, please prepare another PC which has the internet environment.

3 The procedure in the case that the PC where NIS-Elements L is installed is connected to the Internet


In this case, the Online Activation can be executed.

3.1 Online activation

- 1) Please start NIS-Elements L. If the license has not been activated, the following dialog appears. Please click "START" button and move to the next dialog, where Activation ID can be input.

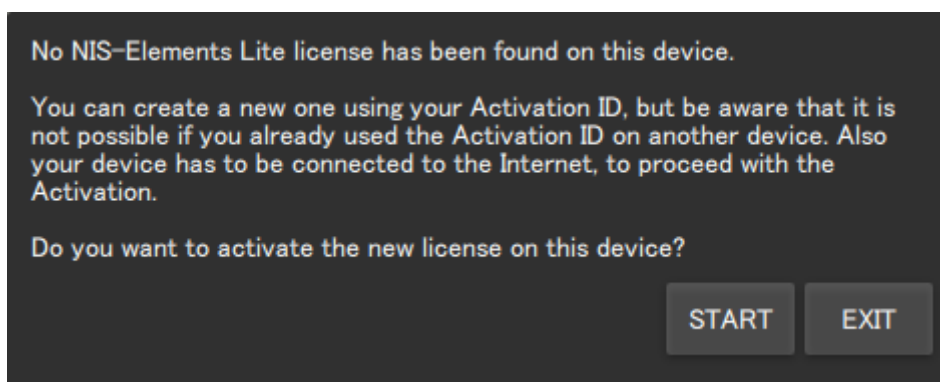


- 2) Please input the activation ID, which is written on the license card, at the following dialog and click "Activate" button. Then, license activation starts.

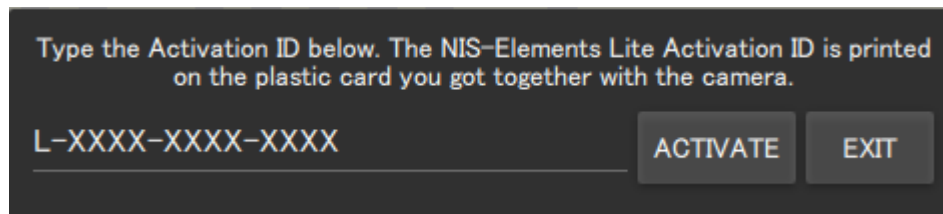


[Note] Please operate this window with Tablet mode in case that hardware keyboard is not connected.

- 3) If activation has been done successfully, the following success dialog is shown. This means license activation is completed. NIS-Elements can be operated after clicking "START" button.

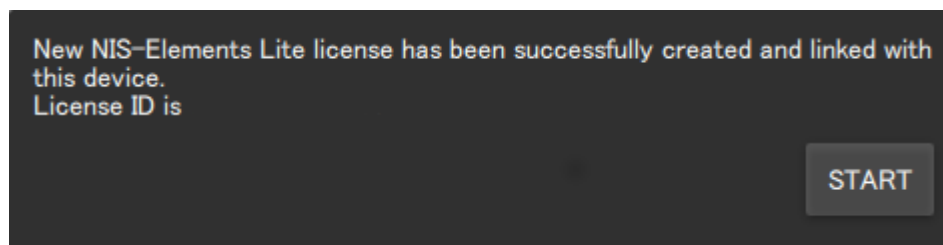


3.2 In the case that online activation is failed

Even if the PC is connected to the Internet, online activation might be failed for some reasons (e.g. not supported proxy server is used). If activation fails, the following error dialog appears. In this case, please follow the below procedure to activate the license.

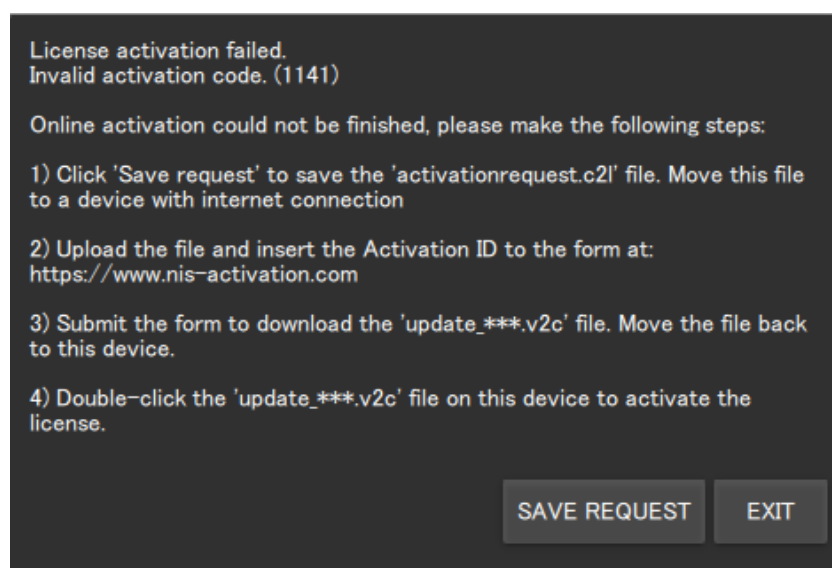

- 1) Please start licmgr.exe, which exists the NIS-Elements L install folder (Default: C:\Program Files\NIS-Elements L). The following window is displayed.

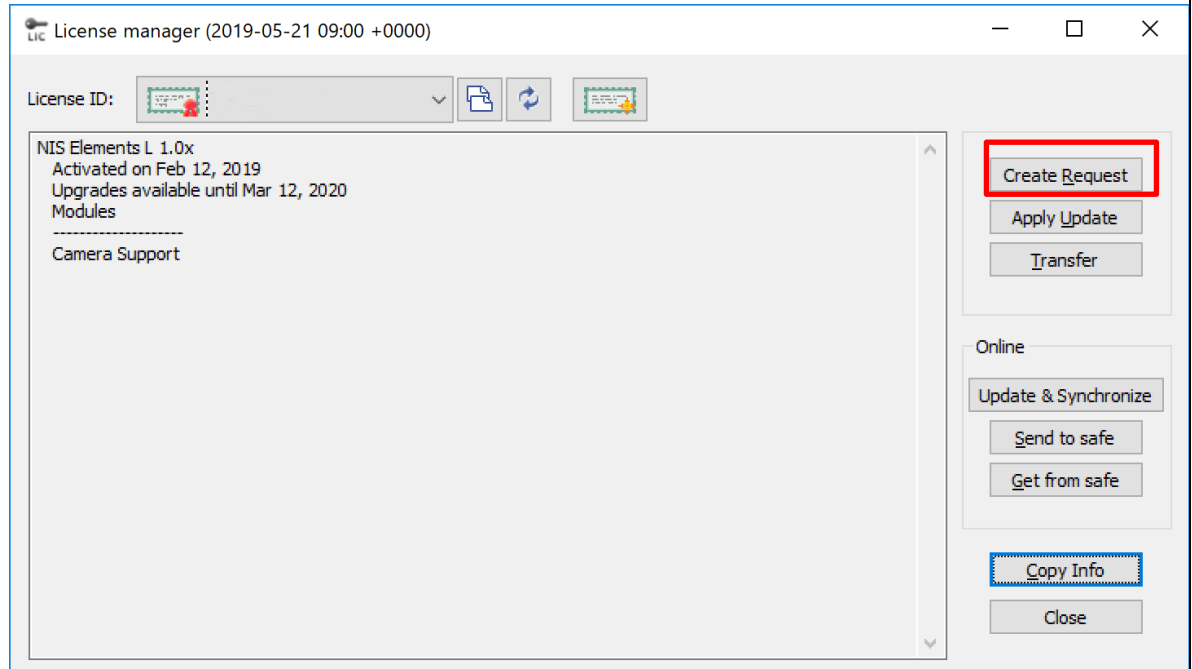


- 2) Please click "Create Request" button and save the request file (C2L file) to any folder. Also, it can be created by "Save Request" button on the error dialog of online activation.

- 3) Start a web browser and access to the following web site.

NIS-Elements L activation site :

<https://www.nis-activation.com/>

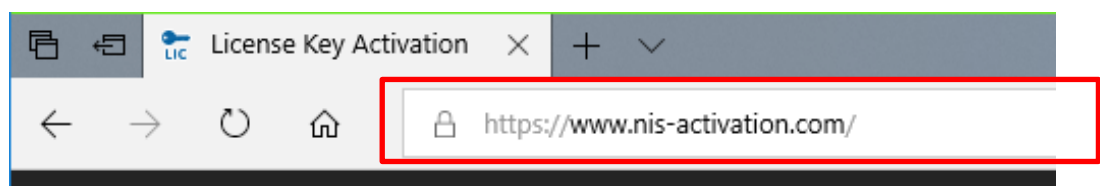

- 4) The following window is shown. Please select the C2L file, created at the step 1, at the "C2L file" and input the Activation ID, written on the license card. Then, click "Create V2C File" and save the activation file, V2C file, to any folder.


License Key Activation

C2L file: 参照...

Activation ID:

Create V2C File

- 5) Please start licmgr.exe again and click "Apply Update" button. Then, select the activation file created at 4) step. After that, the activation is going to be completed. Also, the activation can be completed by double-clicking the v2c file.

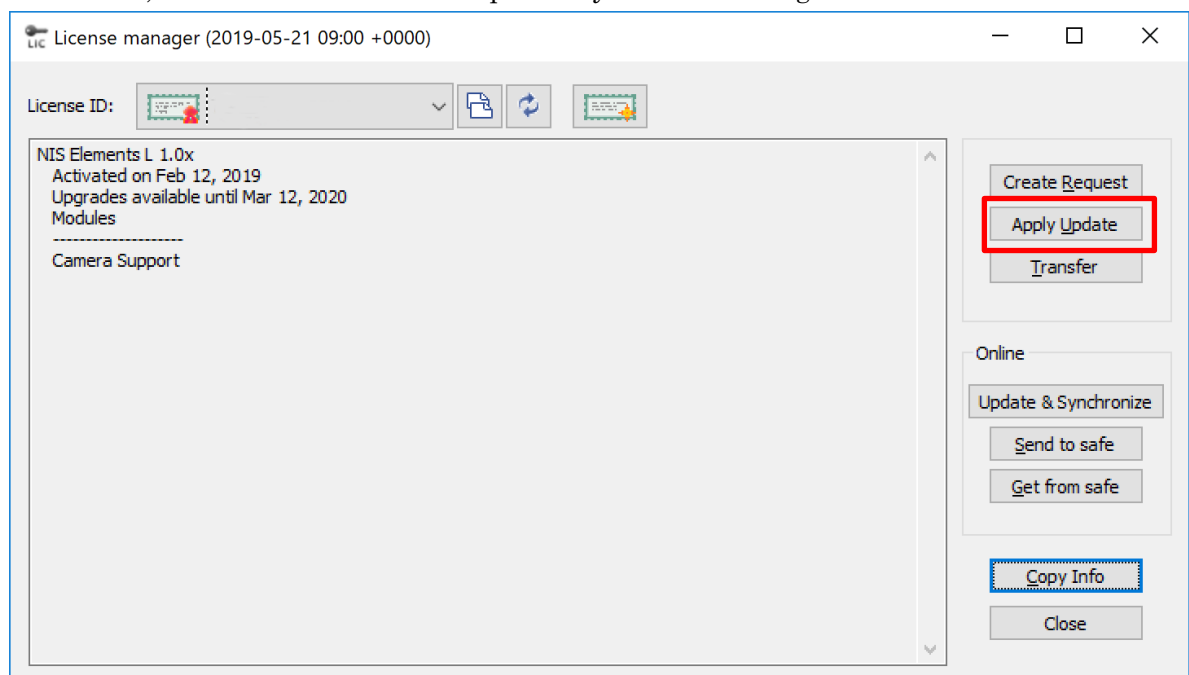


4 In the case that the PC where NIS-Elements L is installed is not connected to the Internet

In this case, another PC which is connected to the Internet is necessary for activating the license. Please follow the below procedure.

4.1 Creating the request file (C2L)

Please perform this step on the PC where NIS-Elements is installed. Start licmgr.exe in the NIS-Elements L install folder (Default: C:\Program Files\NIS-Elements L). Then, click "Create Request" button and save the request file (C2L) to any folder.

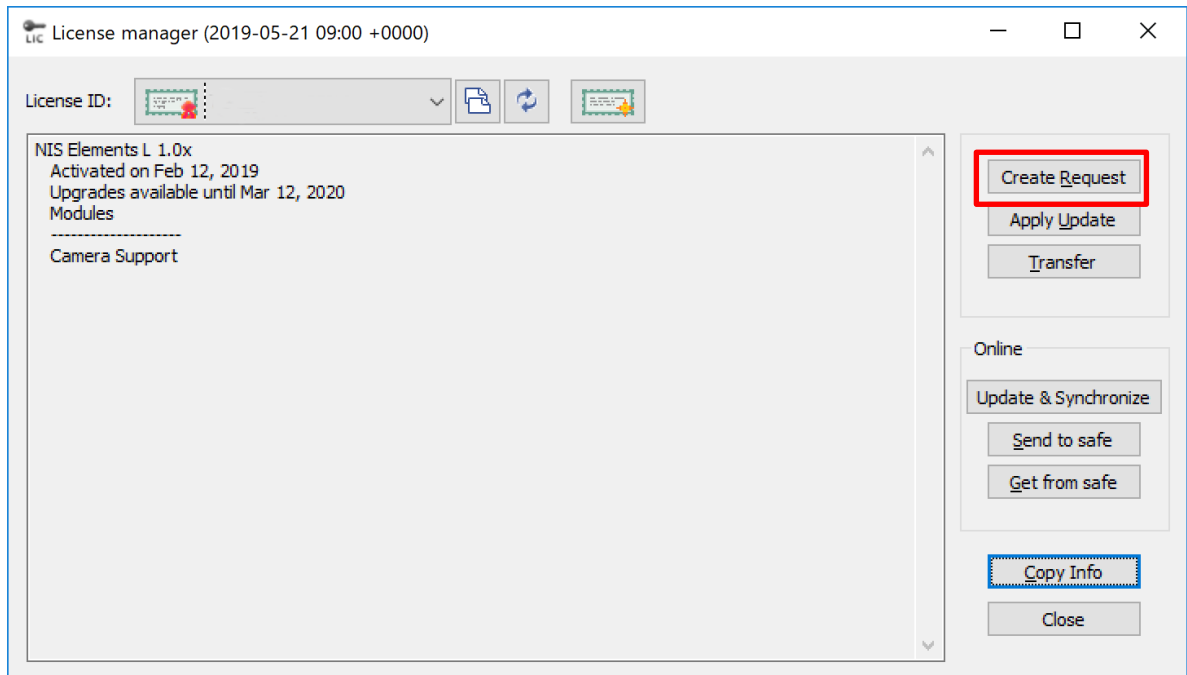


4.2 Creating the activation file (V2L)

Please perform this step on the PC connected to the Internet. First, move the request file (C2L), created at the previous step, to the PC connected to the Internet using USB flash, LAN or etc. Then, start a web browser and access to the following web site.

NIS-Elements L activation site :

<https://www.nis-activation.com/>

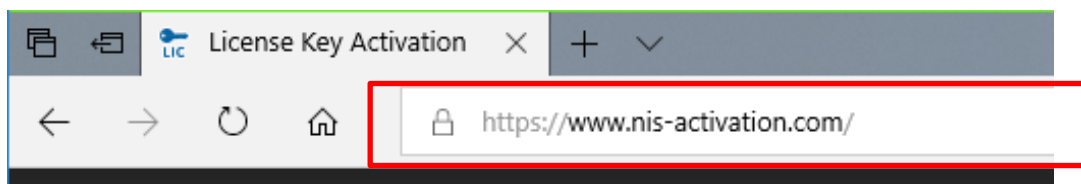

The following window is shown. Please select the C2L file, created at the step 4.1, at the "C2L file" and input the Activation ID, written on the license card. Then, click "Create V2C File" and save the activation file, V2C file, to any folder.


License Key Activation

C2L file:

Activation ID:

4.3 Activating the license

Please perform this step on the PC where NIS-elements L is installed. First, move the V2C file, created at the previous step, to the NIS-Elements L PC using USB flash or LAN. Then, start the licmgr.exe again on the NIS-Elements L PC. Click "Apply update" and load the activation file (V2C), created at the previous step. After that, the activation is going to be completed.

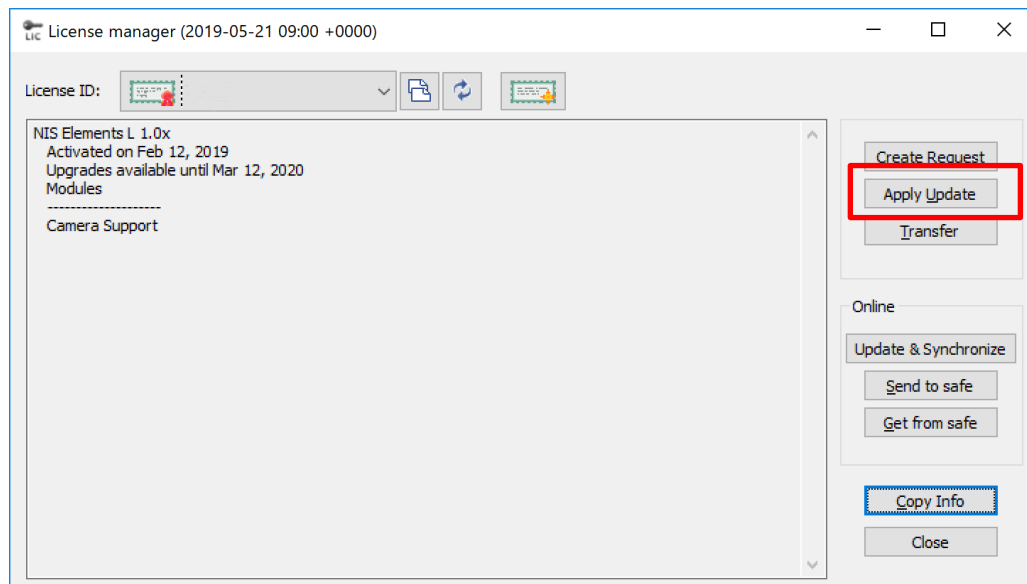

5 [Important] Caution

In case that the PC where license activation is done has to be initialized, it is

necessary for the license to evacuate before it.

The Procedure is described at License Transfer Manual.

End